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Aseptic Heat Exchangers

Calculated and designed individually –
perfectly matching your process

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Operating principle:

The construction of aseptic heat exchangers guarantees the greatest possible separation between the primary and secondary media.

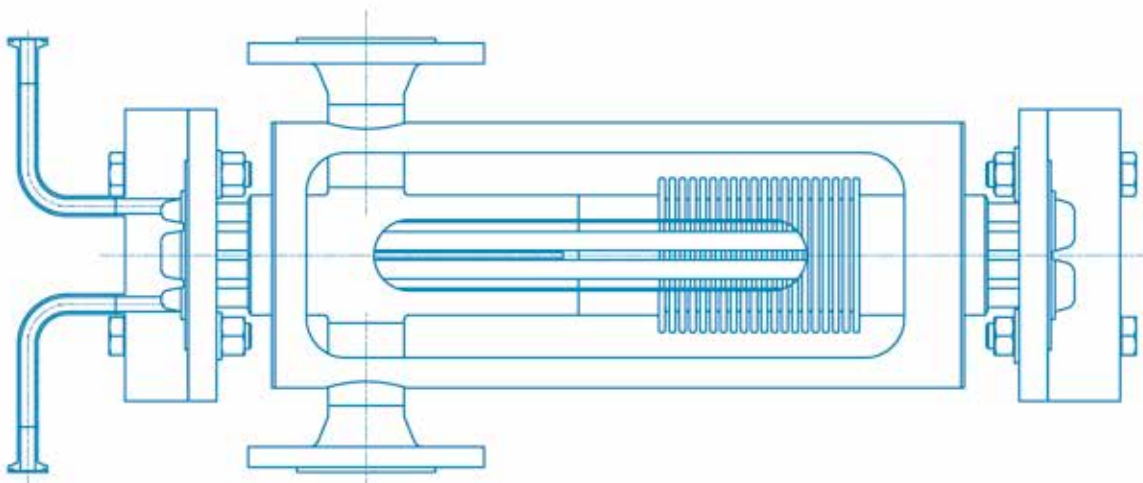
Therefore, the heat exchangers are equipped with a double tubesheet, thus preventing any possibility of mixing and contaminating the processed products with the primary media.

The sterile primary media passes through seamless primary tube via the header and is cooled or heated by the secondary media.

Construction:

NEUMO aseptic heat exchangers are always custom and process specific, thus satisfying the most challenging demands in compliance with your process parameters.

This flexible design reaches highest capacity through adjustment of tube lengths and numbers even if the flow rate is low. To avoid stress cracking caused by differential expansion rates, the outer housing of the heating/cooling chamber is fitted with a sophisticated and highly flexible stainless steel bellow section, guaranteeing operational safety and long life.

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
Features:

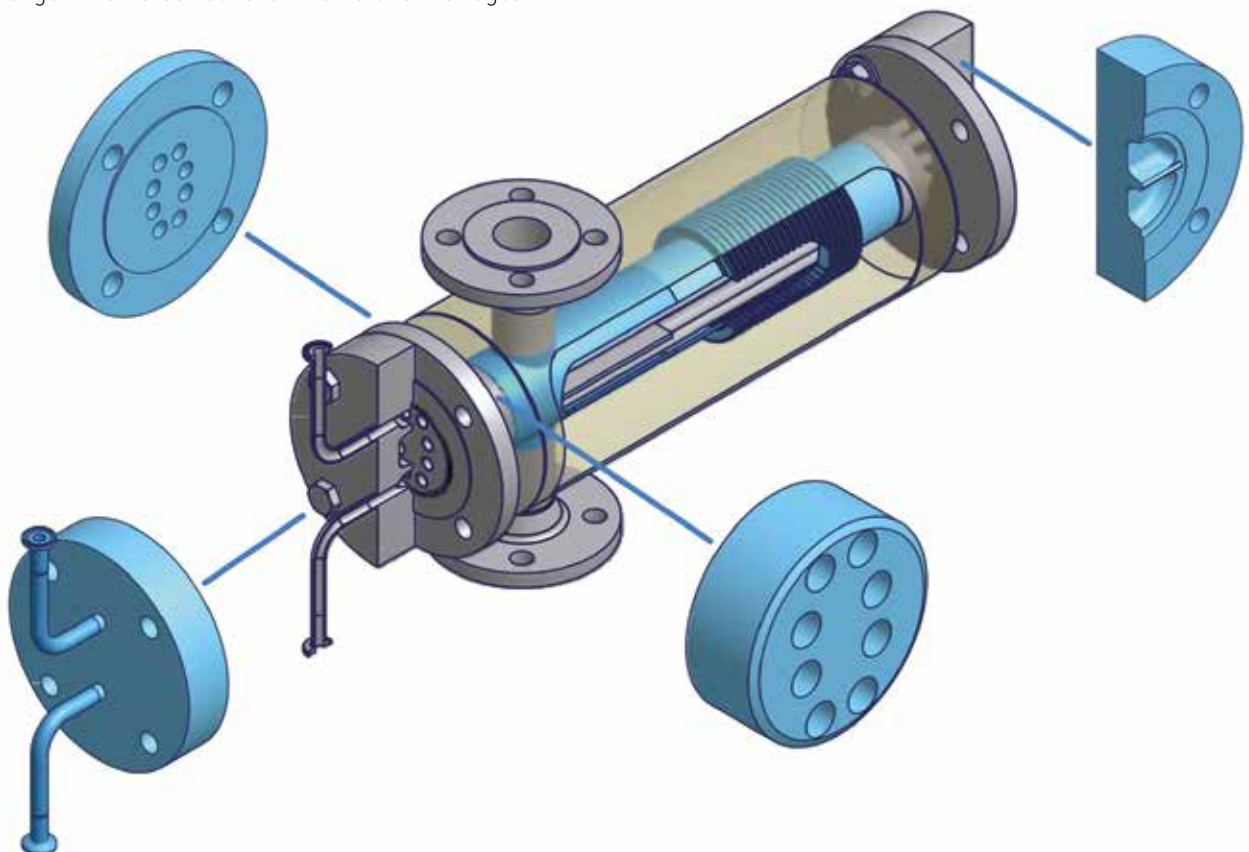- strict separation of processing and cooling / heating media (primary and secondary chamber)
- high pressure and thermal shock resistance due to bellow expansion compensation
- less dead space
- self-draining if correctly mounted
- optimal cleanability / sterilisability (CIP / SIP)
- defined internal surfaces up to $Ra < 0,4 \mu m$ as well as optional electropolishing on the primary side
- quick assembly and disassembly through modular production
- longstanding experience regarding efficient and customised performance design
- dual- or quad-flow construction where space is limited
- horizontal or vertical installation
- qualified sealing material for pharmaceutical areas
- **sealing with NEUMO ConnectS® (elastomer-free!), NEUMO BioConnect®, DIN11864 or according to customer specifications**
- standard material primary circuit: 1.4435 / 1.4404
- special design in nickel-base alloys (such as 2.4602 / 2.4610 / 2.4605)
- insulation available in special materials

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E-Mail: f.titzmann@neumo.de

www.neumo.de

The **NEUMO Ehrenberg Group** is a privately owned and managed multi-national corporate group with more than 2,150 employees.

Founded in 1947, NEUMO, the group's parent company, now produces, manufactures and supplies fittings, equipment and tubes made from stainless steel as well as corrosion-resistant special alloys for the pharmaceutical, biotechnology and chemical industries.

NEUMO products provide a significant contribution to the safety and integrity of production processes.

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NEUMO GmbH + Co. KG